

## C54 PFAS IN SOIL

### 1.0 Sample Reception

- 1.1 All breakages and shortages must be reported within 24 hours of sample receipt.
- 1.2 Store samples at  $4\pm 2^{\circ}\text{C}$  upon receipt.
- 1.3 Check that all the parameters for which you are registered are correctly identified in the PTC portal.
- 1.4 Inquiries regarding samples and their shipment may be directed to:

PT Non-conformances  
Phenova  
Tel: (303) 940-0033  
Email: [AndreaLg@phenova.com](mailto:AndreaLg@phenova.com)

cc: PT Canada Program Officer  
email: [programofficer@PTcanada.org](mailto:programofficer@PTcanada.org)  
cc: Nadine Lewis, PTC Executive Director  
email: [nlewis@PTcanada.org](mailto:nlewis@PTcanada.org)

Inquiries should be made by email. When reporting damage upon receipt, please provide a picture of the damaged samples. Please include your PT Canada laboratory number on all correspondence.

### 2.0 Sample Analysis

- 2.1 Refer to the PTC Catalogue for approximate sample concentrations.
- 2.2 Provided with >1mL spiking solution in methanol in a 2mL amber ampoule; and 30g blank soil in sealed 20mL ampoule.
- 2.3 Assume zero percent moisture for the soil for calculations.
- 2.4 Transfer 10g of the blank soil into your extraction apparatus. Use a gas tight syringe to transfer 50 $\mu\text{L}$  of the spiking solution onto the 10g blank soil. Extract and proceed with testing using the routine analytical method identified in your PT Canada application.
- 2.5 Each sample contains random selection percentage of analytes with at least 60% of the PFAS analytes in each sample. All analytes are covered per PT Study. Please see the list of PFS analytes below.

### 3.0 Reporting Results

- 3.1 Results must be reported by midnight of the study deadline in the PTC portal.
- 3.2 Report RDL (optional) if you want RDL accounted for in z scores.

### 4.0 Safety

- 4.1 The PT samples are designed for use by laboratory professionals familiar with environmental samples and potentially hazardous materials.

## **PFAS, EPA 1633 Analyte List (40 Analytes)**

| <b>Analyte Description</b>             | <b>CAS Number</b> |
|----------------------------------------|-------------------|
| Perfluorobutanoic acid (PFBA)          | 375-22-4          |
| Perfluoropentanoic acid (PFPeA)        | 2706-90-3         |
| Perfluorohexanoic acid (PFHxA)         | 307-24-4          |
| Perfluoroheptanoic acid (PFHpA)        | 375-85-9          |
| Perfluorooctanoic acid (PFOA)          | 335-67-1          |
| Perfluorononanoic acid (PFNA)          | 375-95-1          |
| Perfluorodecanoic acid (PFDA)          | 335-76-2          |
| Perfluoroundecanoic acid (PFUnA)       | 2058-94-8         |
| Perfluorododecanoic acid (PFDoA)       | 307-55-1          |
| Perfluorotridecanoic acid (PFTriA)     | 72629-94-8        |
| Perfluorotetradecanoic acid (PFTeA)    | 376-06-7          |
| Perfluorobutanesulfonic acid (PFBS)    | 375-73-5          |
| Perfluoropentanesulfonic acid (PFPeS)  | 2706-91-4         |
| Perfluorohexanesulfonic acid (PFHxS)   | 355-46-4          |
| Perfluoroheptanesulfonic Acid (PFHpS)  | 375-92-8          |
| Perfluorooctanesulfonic acid (PFOS)    | 1763-23-1         |
| Perfluorononanesulfonic acid (PFNS)    | 68259-12-1        |
| Perfluorodecanesulfonic acid (PFDS)    | 335-77-3          |
| Perfluorododecanesulfonic acid (PFDoS) | 79780-39-5        |
| Perfluorooctanesulfonamide (FOSA)      | 754-91-6          |
| NMeFOSAA                               | 2355-31-9         |

|                |             |
|----------------|-------------|
| NEtFOSAA       | 2991-50-6   |
| 4:2 FTS        | 757124-72-4 |
| 6:2 FTS        | 27619-97-2  |
| 8:2 FTS        | 39108-34-4  |
| NEtFOSA        | 4151-50-2   |
| NMeFOSA        | 31506-32-8  |
| NMeFOSE        | 24448-09-7  |
| NEtFOSE        | 1691-99-2   |
| 9Cl-PF3ONS     | 756426-58-1 |
| HFPO-DA (GenX) | 13252-13-6  |
| 11Cl-PF3OUdS   | 763051-92-9 |
| ADONA          | 919005-14-4 |
| 3:3 FTCA       | 356-02-5    |
| 5:3 FTCA       | 914637-49-3 |
| 7:3 FTCA       | 812-70-4    |
| NFDHA          | 151772-58-6 |
| PFMBA          | 863090-89-5 |
| PFMPA          | 377-73-1    |
| PFEESA         | 113507-82-7 |